



Nutrient Analysis, Nitrate, Nitrite, Silicate, Phosphate and Ammonium

SUMMARY: The phytoplankton macro nutrients nitrate, nitrite, silicate, and phosphate in seawater are analyzed using colorimetric assays. Ammonium concentrations are determined using a fluorometric assay.

1. Principle

Nutrient analysis is performed on a QuAAtro continuous flow autoanalyzer (SEAL Analytical). A sample of seawater enters a reagent stream within a manifold on the analyzer where it undergoes a series of reactions that ultimately produce a colored compound. These compounds absorb light at a specific wavelength. A monochromatic beam of light is passed through the sample and the absorbance is measured. The machine is calibrated with a series of known standards and a standard curve is produced. The intensity of the color produced by the unknown sample is proportional to the concentration of the analyte present. The product of the ammonia method is a fluorescent species; however the same basic principle applies, where the intensity of the fluorescence is directly related to concentration. The methods for silicate and total oxidized nitrogen (TON) are modified versions of those described by used Armstrong et al. (1967) and Gordon et al. (1992). The phosphate determination employs a modification of the method described by Bernhardt and Wilhelms (1967), and ammonia is analyzed based on the Kerouel and Aminot (1997) fluorometric method.

2. Method Description

Silicate (SiO₂)

Silicate is analyzed using a modified technique of Armstrong (1967). An acidic solution of ammonium molybdate is added to a seawater sample to produce silicomolybdic acid which is then reduced to a blue silicomolybdous acid following the addition of ascorbic acid. The amount of blue color produced is proportional to the amount of dissolved reactive silicate in the sample. Oxalic acid is added to inhibit PO₄ color interference. The sample is passed through a 10mm flow cell and the absorbance is measured at 820nm.

Nitrate and Nitrite (NO₃ and NO₂)

A modification of the Armstrong (1967) procedure is used for the analysis of nitrate plus nitrite, or total oxidized nitrogen (TON). For this analysis, the seawater sample is passed through a cadmium reduction coil where nitrate is reduced to nitrite. The efficiency of this reduction is determined by running two equimolar solutions, one containing only nitrate and one containing only nitrite, through the coil. The percent of nitrate converted to nitrite yields the coil efficiency, a factor used to calculate the nitrate concentration. Sulfanilamide is introduced to the sample stream followed by N-(1-naphthyl) ethylenediamine dihydrochloride which complexes with nitrite to form a red azo dye. The stream is then passed through a 10 mm flowcell and the absorbance measured at 520nm.

The same method is employed for nitrite analysis, except the cadmium column is not present. Consequently, only nitrite already present in the sample is measured.

The nitrate concentration is calculated on a "virtual channel" by the following equation, using the coil

efficiency, TON, and NO₂:

$$NO_3 = \frac{\left[TON * \frac{(A * Recovery + B)}{A + B} \right] - NO_2}{Recovery}$$

NO₃=Nitrate (in sample)

NO₂= Nitrite (in sample)

TON=Total oxidized nitrogen (NO₃+NO₂) in sample

A= NO₃ concentration in mixed calibrant

B= NO₂ concentration in mixed calibrant

Recovery = Coil efficiency (expressed as a decimal)

Phosphate (PO₄)

Phosphate is analyzed using a modification of the Bernhardt and Wilhelms (1967) technique. Similar to the silicate method, an acidic solution of ammonium molybdate is added to the sample to produce phosphomolybdic acid that is subsequently reduced to blue phosphomolybdous acid following the addition of ascorbic acid. Color intensity is directly related to the concentration of dissolved phosphate in the sample. The reaction product is then passed through a 10mm flow cell and the absorbance measured at 880nm.

Ammonia (NH₃)

Ammonia is measured fluorometrically using a modification of the method described by Kerouel and Aminot (1997). In the presence of a borate buffer, samples are reacted with o-phthalaldehyde (OPA) to form a fluorescent complex that is excited at 370nm and emits at 460nm. The reaction takes place at 75°C. Sodium sulfite is added to the working reagent to reduce sensitivity to dissolved amino acids.

3. Method Notes

Ammonium is a difficult parameter to measure accurately due to its insidious nature and problems with contamination. Care must be taken during sampling to insure there is no contamination (e.g. touching the inside of the tube or the cap with fingers, smoking near rosette).

Standards are made up in artificial seawater comprised of 34g NaCl and 0.5g NaHCO₃ per liter of ultrapure water. Batches are generally only used for 2-3 days and then discarded to minimize the potential for ammonia contamination.

4. Water Sampling

Nutrient samples are drawn into 50 ml polypropylene screw-capped centrifuge tubes.

The tubes and caps are cleaned with 10% HCl and rinsed 3 times with sample before filling.

Samples are analyzed within 2-16 hours after sample collection, allowing sufficient time for all samples to reach room temperature. The centrifuge tubes fit directly onto the sampler.

5. Calculations

All data is reported in micro-moles/liter. The main calculations for concentration on the Bran and Luebbe AA3 are run through the required AACE software interface. These calculations still follow the principle of other instruments, where:

$$[X] \text{ micro moles/liter} = (\text{Absorbance-blank}) \times F1 \text{ (Response Factor)}$$

Values are corrected for drift based on changes in beginning and end standards, ultra pure water and the relative position of samples in the run. Corrections for linearity are performed, if necessary, based on a set of absorbences and concentrations; deviations from Beer's law can be plotted to reveal a polynomial function that can be applied to correct sample values accordingly. Improvements in optics in the AA3 & QuAAtro instruments have resulted in marked improvement in linearity and reduction of blank values for nitrate and silicate and phosphate.

The nature of ammonium poses a particular problem in relation to reliable blank values. Upon expert recommendation, the true baseline blank is based on the value obtained from deep seawater samples (500m) where ammonium is thought to be zero. Carefully monitoring these values in relation to machine function enables them to be used as ammonium blanks and used to correct all depths.

6. Quality Control

A sample of reference material for nutrients in seawater (RMNS), produced by KANSO technos (www.kanso.co.jp) is included in every run and those data are used to adjust values for nitrate, nitrite, phosphate, and silicate if appropriate.

An aliquot from a large volume of stable deep seawater is run with each set of samples as an additional check. The stability of the deep seawater check is aided by the addition of mercuric chloride as a poison.

The efficiency of the cadmium column used for nitrate reduction is monitored throughout the cruise and usually ranges from 98.0-100.0%.

NO_3 , PO_4 , NO_2 , and NH_4 are reported to two decimal places and SiO_2 to one.

Accuracy is based on the quality of the standards; the levels in micro moles/liter (μM) are:

$$\text{NO}_3 = 0.05$$

$$\text{PO}_4 = 0.004$$

$$\text{SiO}_2 = 2-4$$

$$\text{NO}_2 = 0.05$$

$$\text{NH}_3 = 0.03$$

The precision of the instrument for NO_3 , PO_4 , and NH_4 is 0.01 μM and 1.0 μM for silicate

and 0.01 μM for NO_2 .

The detection limits in micro moles/liter for the instrumentation are:

$$\text{NO}_3 + \text{NO}_2 = 0.02$$

$$\text{PO}_4 = 0.02$$

$$\text{SiO}_2 = 0.5$$

$$\text{NO}_2 = 0.02$$

$$\text{NH}_3 = 0.04$$

7. Equipment/Supplies

Seal Analytical continuous-flow QuAatro (run by IOD since CalCOFI 1203SH; AutoAnalyzer 3 (AA3) run by ODF on cruises prior to 1203SH. Distributed by Bran and Luebbe, <http://www.seal-analytical.com/>

50 ml centrifuge tubes and test tube racks, 6 sets color coded and numbered

Milli-Q purified water system or equivalent polished water source

Sundry laboratory glassware

8. References

Armstrong, F.A.J., Stearns, C.A., and Strickland, J.D.H., (1967). "The measurement of upwelling and subsequent biological processes by means of the Technicon Autoanalyzer and associated equipment," Deep-Sea Research, 14, pp.381-389.

Atlas, E.L., Hager, S.W., Gordon, L.I., and Park, P.K., (1971). "A Practical Manual for Use of the Technicon AutoAnalyzer in Seawater Nutrient Analyses Revised," Technical Report 215, Reference 71-22, p.49, Oregon State University, Department of Oceanography.

Bernhardt, H., and Wilhelms, A., (1967). "The continuous determination of low level iron, soluble phosphate and total phosphate with the AutoAnalyzer," Technicon Symposia, 1, pp.385-389 .

Gordon, L.I., Jennings, J.C., Ross, A.A., Krest, J.M., (1992). "A suggested Protocol for Continuous Flow Automated Analysis of Seawater Nutrients in the WOCE Hydrographic Program and the Joint Global Ocean Fluxes Study," Grp. Tech Rpt 92-1, OSU College of Oceanography Descr. Chem Oc. .

Hager, S.W., Atlas, E.L., Gordon L.I., Mantyla, A.W., and Park, P.K., (1972). " A comparison at sea of manual and autoanalyzer analyses of phosphate, nitrate, and silicate ," Limnology and Oceanography, 17, pp.931-937.

Keuroul, R., Aminot, A., (1997) "Fluorometric determination of ammonia in sea and estuarine waters by direct segmented flow analysis," Marine Chemistry Vol. 57, no. 3-4, pp. 265-275.

Last Updated on Wednesday, 28 August 2013 10:44